# Mycoplasma Removal Agent

## (Cell Culture Grade)

### Ordering info

TBR0332. Mycoplasma Removal Agent 50x, 100 mL

### Description

Mycoplasma is the predominant contaminant in cell cultures, posing challenges in detecting pollutants due to interference with experimental results. Cell culture mycoplasma contamination is estimated to occur in approximately 30.3% to 50.5% of cases.

**Mycoplasma Removal Agent, 50x** is an antibiotic solution based on a derivative of the quinoline. It is a highly effective against multiple mycoplasma species. Quinoline compound eliminates mycoplasma infection by inhibiting mycoplasma DNA gyrase.

### Storage

Store at -20°C. Stable for 2 years.

The product is shipped on blue ice.

### Features

- Ready to use.
- Low cytotoxicity for eucaryotic cell.
- Endotoxin  $\leq 10$  EU/ mL.

### Applications

- Elimination of multiple mycoplasma species.
- Prevention of contamination of cultures at 0.1  $\mu\text{g}/\text{mL}$ .

Also available:

[Resazurin Viability Assay Kit \(TBK0506, TBK0507\)](#)

[G-418 Sulfate Solution 50 mg/mL \(TBR0334\)](#)

[Trypan Blue Cell Viability Indicator \(TBB0402\)](#)


[Trypsin/ EDTA \(TBZ0342, TBZ0344\)](#)

[PBS 1x, pH=7.4 \(TBB0360\)](#)

## PROTOCOL

*Mycoplasma removal includes 3 or more passages. It is important to monitor weekly for mycoplasma presence. Contamination should be removed after 2-3 weeks. Once the cells are mycoplasma free, transfer them to a growth media without Mycoplasma Removal Agent and confirm that mycoplasma contamination has been eliminated.*

1. Carefully, aspirate all the culture media off.
2. Wash the cells with prewarmed PBS 1x.
3. Detach cells using Trypsin/ EDTA (TBZ0342, TBZ0344) or Accutase (TBZ0340).
4. Count the cells.
5. Seed the cells in fresh medium supplemented with Mycoplasma Removal Agent (1 mL per 49 mL of growth media).
6. Incubate the cells in appropriate conditions for 2-3 days.

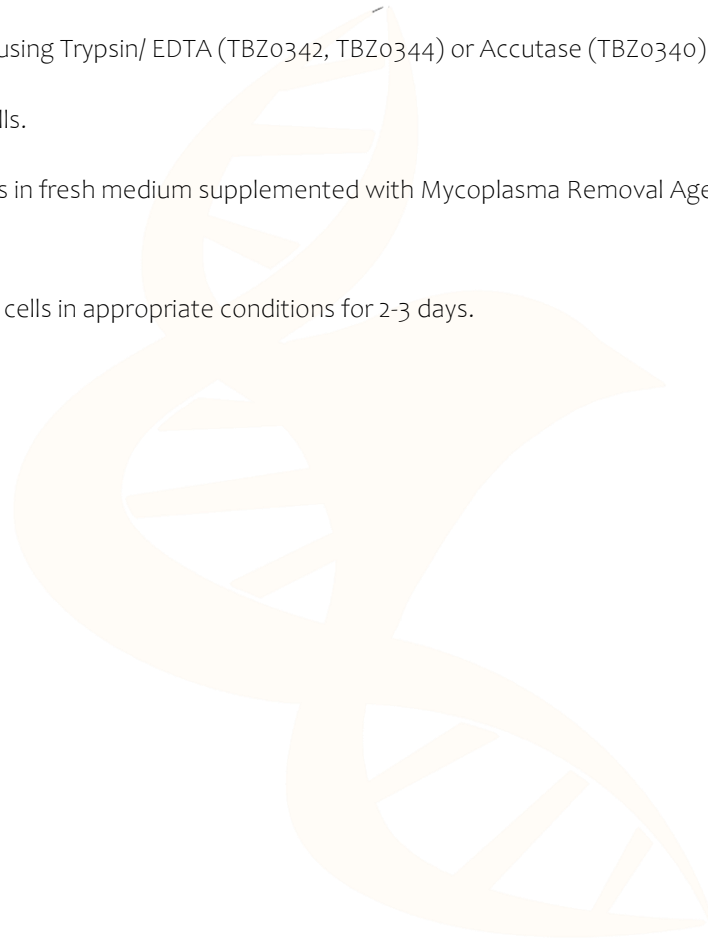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Resazurin Viability Assay Kit (TBK0506, TBK0507)

G-418 Sulfate Solution 50 mg/mL (TBR0334)

Trypan Blue Cell Viability Indicator (TBB0402)

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