Carbenicillin

(Disodium Salt)

Ordering info

TBR0119, Carbenicillin 1 g

Description

Carbenicillin Disodium Salt is a semi-synthetic penicillin antibiotic with a relatively greater spectrum of activity against Gram-negative bacteria. Carbenicillin binds and inhibits the enzymes involved in the synthesis of bacterial cell wall. Since carbenicillin is less sensitive to β -lactamases than ampicillin, it is effective against ampicillin-resistant bacteria. It has superior stability at low pH and acts as a selection agent.

CAS: 4800-94-6

Formula: $C_{17}H_{16}N_2Na_2O_6S$

MW: 422.36 g/mol

Synonym: α -Carboxybenzylpenicillin

Features

- Highest purity.
- Hygroscopic white powder.
- Solubility in water: 50-100 mg/mL.
- Higher stability than ampicillin.

Storage

Store at 2-8°C.


Following reconstitution, aliquot and freeze at -20°C.

Quality Control

Purity is tested by HPLC (>90 %).

Applications

- Used to select for ampicillin resistance in mutated and transformed cells (preferred over ampicillin due to its increased stability).
- Commonly used for the elimination of *Agrobacterium* species after inoculation.

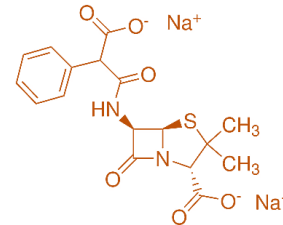PROCEDURE TO PREPARE CARBENICILLIN STOCK SOLUTION (50 mg/ mL)

1. Dissolve **1g Carbenicillin** (TBR0119) in water. Add **water up 20 mL**.
2. Prewash a 0.22 – 0.45 μ m filter with 20 mL of water.
3. Put the washed filter over a new sterile tube. Then, filter Carbenicillin Stock Solution.
4. Make aliquots in 1.5 mL tubes and store at -20°C.

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